

Supercritical Fluid Extraction of Residual Pesticides in Coffee Bean

Introduction

In Japan on the 29th of May 2006 the Ministry of Health, Labor and Welfare (MHLW) promulgated the Positive List System for residual pesticides, food additives, and veterinary medicines remaining in foods, following the revision of the Food Hygiene Law. In this list approximately 800 kinds of those agricultural chemicals were registered. This system is to prohibit the distribution of foods that contain more than 0.01 ppm of each chemical.

The extraction of residual pesticides in foods has been performed by the solvent extraction method. This method, however, takes about 4 - 5 hours for each extraction, and requires a huge amount of organic solvent. In recent years, supercritical fluid extraction (SFE) using supercritical carbon dioxide has attracted much attention as an alternative method to the solvent extraction method.

We have developed a fully automated residual pesticide extraction system, and applied this system to the analysis of a coffee bean sample. Extracted components were analyzed by GC-MS/MS.

Experimental

The newly developed fully automated residual pesticide extraction system was used throughout the experiment. The schematic diagram of this system is shown in Figure 1.

A sample of coffee beans was selected for analysis. Sixty-eight kinds of pesticides were added to the coffee beans to be a concentration of 0.1 ppm for each pesticide except captan, 1 ppm and acetamiprid, 0.5 ppm. Three grams of the coffee beans was loaded in each extraction vessel; SFE was applied at an extraction pressure of 15 MPa, at an extraction temperature of 40^o C, for an extraction time of 30 min; the extracted components were adsorbed on a trap column; the trapped components were eluted with acetonitrile; the acetonitrile solution was evaporated to dryness with nitrogen gas; and the residue was dissolved in 3 mL of acetone containing 0.05% of PEG200 and PEG400. A portion of this solution was injected onto the GC.

Results and Discussion

Chromatograms of the standard mixture (upper), the sample added with the standard (middle), and the blank (lower) are shown in Figure 2.

As shown in Table 1, among 68 components of the pesticides, 58 components exhibited more than 70% recovery, and 66 components more than 50% recovery. The recovery of acetamiprid and penycuron was as low as 42% and 14%, respectively.

Acetamiprid, owing to its high hydrophilicity, indicated low solubility in supercritical carbon dioxide, resulting in a poor recovery in SFE. Low recovery of penycuron seemed to be ascribed to the sample matrix.

References

- 1) Ministry of Health, Labor and Welfare Official Gazette No. 498
- 2) Ministry of Health, Labor and Welfare Official Gazette No. 497

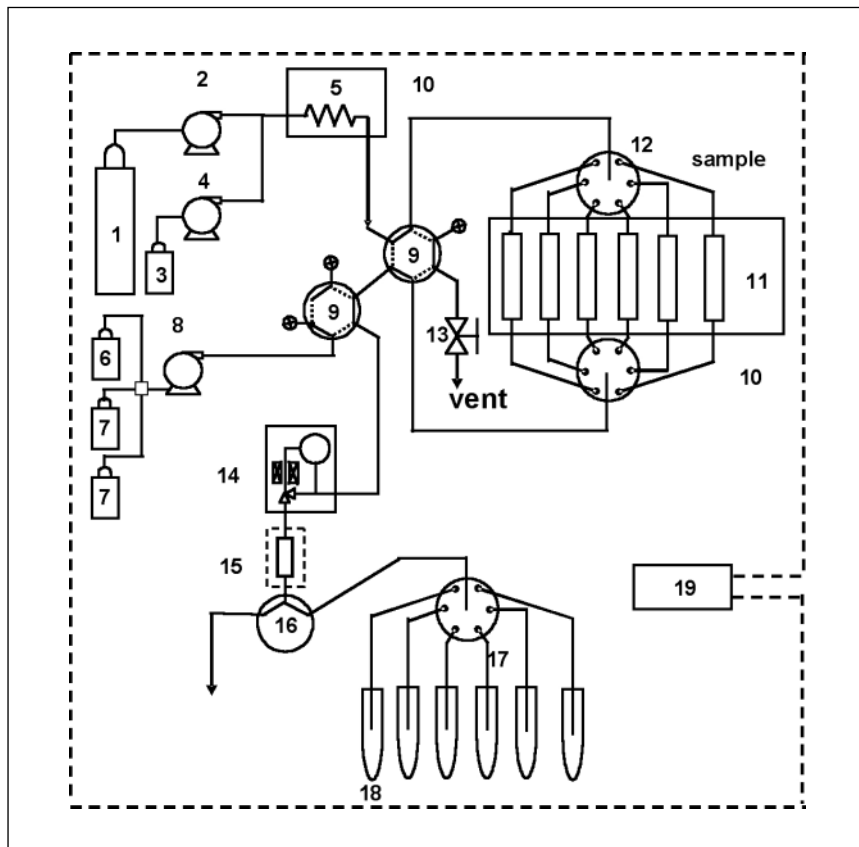


Figure 1 Schematic Diagram of fully automated system for supercritical fluid extraction of residual pesticides

System configuration: 1 = carbon dioxide cylinder, 2 = liquefied carbon dioxide delivery pump, 3 = modifier, 4 = modifier delivery pump, 5 = preheating coil, 6 = solvent for trap elution, 7 = rinse solution for trap column, 8 = solvent delivery pump, 9 = switching valve for flow line, 10 = oven, 11 = extraction vessels, 12 = 6-vessel changer, 13 = release valve, 14 = automatic back pressure regulator, 15 = trap column, 16 = 3-way valve, 17 = 6-way flow line switching valve, 18 = collection tubes, 19 = system controller

Supercritical fluid extraction conditions: extraction tube = 10 mL (10 mm x 127 mm), supercritical fluid = CO₂, back pressure = 15 MPa, extraction time = 30 min, flow rate = 2 mL/min, trap column = ODS (4.6 mm x 50 mm, 30 μm), solvent for trap elution = acetonitrile 2 mL (flow rate = 2 mL/min).

Supercritical CO₂ delivered by pump 2 passes through one of the vessels 11 in which the sample is loaded and then pesticides are extracted. The extracted pesticides are concentrated by the trap column, is eluted by acetonitrile (2 mL) delivered by pump 8, and is collected in one of collection tubes 18. This system is automatically controlled by 19, system controller.

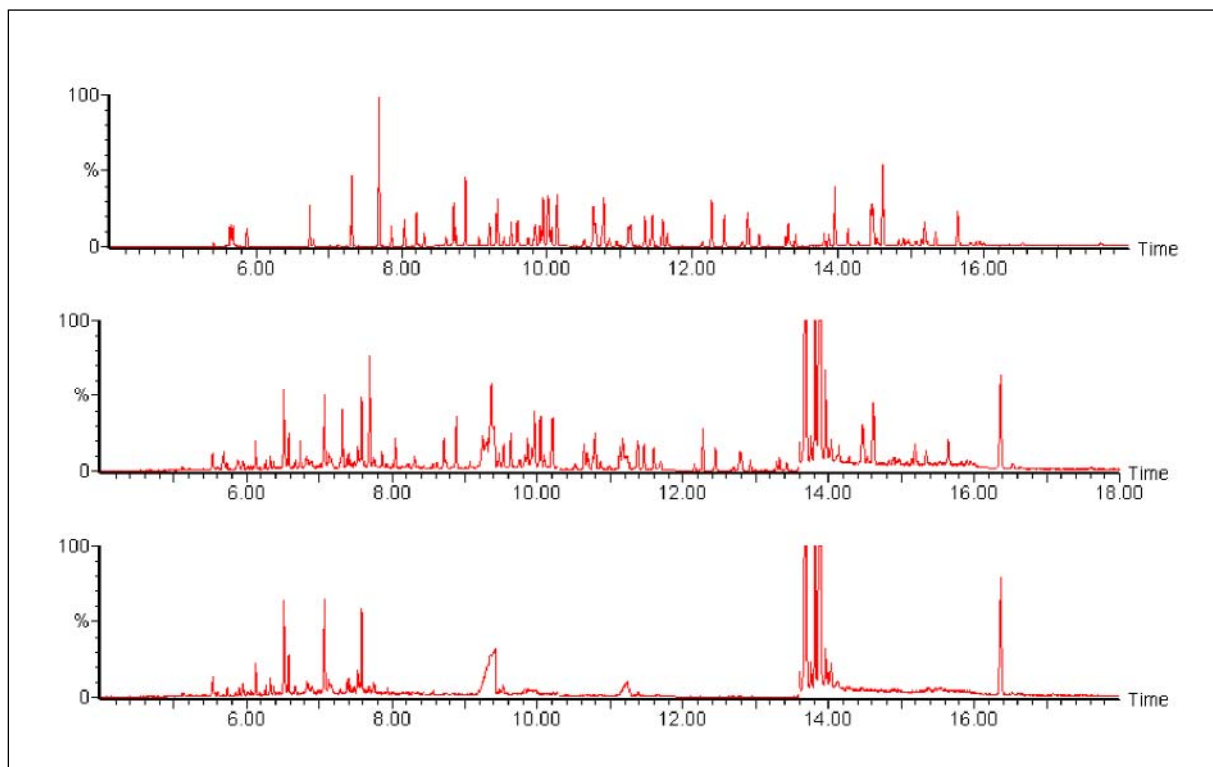


Figure 2 GC chromatograms of wheat flour sample. Upper: standard mixture (68 components), Middle: sample added with standard mixture, and Lower: blank

Measurement conditions: Instrument=Quattro micro GC (Waters micromass), Ionization method = EI, Measurement mode=MRM, SIM, Ionization source temperature=280 °C, Interface temperature=280 °C, GC=6890N(Agilent), Injection method=Splitless, Injection volume=1 μL, Inlet temperature=250° C, Column=DB-5MS(30 m x 0.25 mm), Column temperature=50 °C (0 min)—50 °C (1 min) — 200 °C (7 min) — 250 °C (9 min) — 300 °C (11 min).

Standard mixture solution contains 68 components as below.: 1: Acephate, 2: Acetamiprid, 3: Bendiocarb, 4: Bitertanol, 5: Butylate, 6: Captan, 7: Carbaryl 8: Chinomethionat, 9: Chlorfenvinphos, 10: Chlorpyriphos, 11: Cyfluthrin, 12: Cypermethrin, 13: Deltamethrin, 14: Diazinon, 15: Dichlofluanid, 16: Dichlorvos, 17: Diethofencarb, 18: Dimethylvinphos, 19: EPN, 20: Esprocarb, 21: Ethiofencarb, 22: Ethoprophos, 23: Fenarimol, 24: Fenitrothin, 25: Fenobucarb, 26: Fensulfothion, 27: Fenvalerate, 28: Flucythrinate, 29: Flusilazole, 30: Flutolanil, 31: Fluvalinate, 32: Flutoranil, 32: Imibenconazole, 33: Iprodione, 34: Isofenphos, 35: Isofenphos P=O, 36: Isoprocarb, 37: Lenacil, 38: Malathion, 39: Mefenacet, 40: Mepronil, 41: Methamidophos, 42: Metolachlor, 43: p,p'-DDE, 44: Paclobutrazol, 45: Pencycuron, 46: Pendimethalin, 47: Permethalin, 48: Phenthoate, 49: Phosalone, 50: Pirimifos-methyl, 51: p,p'-DDD, 52: Pretilachlor, 52: Pretilachlor, 53: Propiconazole, 54: Pyraclofos, 55: Pyridaben, 56: Pyridaphenthion, 57: Pyrimidifen, 58: Quinalphos, 59: Tefluthrin, 60: Terbucarb, 62: Thenylchlor,

63: Tolclofos-methyl, 64: Triadimenol, 65: α-BHC, 66: β-BHC, 67: γ-BHC, 68: δ-BHC .

Table 1 The recovery of each pesticide

No	Pesticide	Recovery (%)	No	Pesticide	Recovery (%)	No	Pesticide	Recovery (%)
1	Acephate	95.4	24	Fenitrothion	91.8	47	Permethrin	73.3
2	Acetamiprid	42.2	25	Fenobucarb	77.6	48	Phenthoate	67.6
3	Bendiocarb	84.2	26	Fensulfothion	88.9	49	Phosalone	90.2
4	Bitertanol	81.8	27	Fenvalerate	79.6	50	Pirimifos-methyl	78.7
5	Butylate	62.6	28	Flucythrinate	83.3	51	p,p'-DDD	76.2
6	Captan	54.6	29	Flusilazole	73.5	52	Pretilachlor	75.6
7	Carbaryl	99.1	30	Flutolanil	101.3	53	Propiconazole	73.5
8	Chinomethionat	94.6	31	Fluvalinate	88.7	54	Pyraclufos	76.9
9	Chlorfenvinphos	73.0	32	Imibenconazole	70.3	55	Pyridaben	72.1
10	Chlorpyriphos	80.5	33	Iprodione	92.0	56	Pyridaphenthion	85.0
11	Cyfluthrin	85.6	34	isofenphos	57.7	57	Pyrimidifen	71.0
12	Cypermethrin	78.4	35	Isofenphos P=O	79.5	58	Quinalphos	72.0
13	Deltamethrin	97.3	36	Isoprocarb	70.8	59	Tefluthrin	70.2
14	Diazinon	72.9	37	Lenacil	63.8	60	Terbucarb	102.1
15	Dichlofluanid	85.6	38	Malathion	91.9	61	Terbufos	65.6
16	Dichlorvos	61.6	39	Mefenacet	85.4	62	Thenylchlor	73.3
17	Diethofencarb	80.6	40	Mepronil	75.7	63	Tolclofos-methyl	73.9
18	Dimethylvinphos	71.9	41	Methamidophos	65.4	64	Triadimenol	71.3
19	EPN	71.2	42	Metolachlor	73.5	65	α -BHC	74.5
20	Esprocarb	109.5	43	p,p'-DDE	70.0	66	β -BHC	73.3
21	Ethiofencarb	104.4	44	Paclobutrazol	112.7	67	γ -BHC	70.9
22	Ethoprophos	74.4	45	Pencycuron	14.3	68	δ -BHC	78.1
23	Fenarimol	74.3	46	Pendimethalin	74.4			