
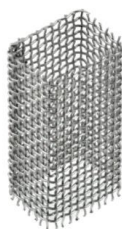


Special Test Accessories

Sometimes drug monographs require modified equipment that is different to the standard Paddle or Basket method. We can provide this for the following:

Felodipine This uses a special suspended basket that is fixed to a modified vessel cover. This is available for most baths.

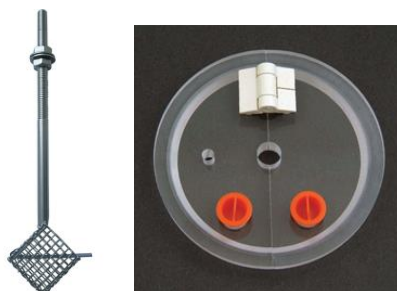

Metformin The 500/1000mg versions of the tablet require special quadrangular inserts inside the standard USP baskets to enable the capsule to remain vertical during the test. These work in any bath.



Alfuzosin Hydrochloride The extended release version uses a special cylindrical sinker. These work in any bath.



Dabigatran The 150mg capsule uses a wider 24.5mm basket. This basket requires a special basket shaft to hold it in place which is available for most dissolution baths. This method was developed by LabHut/QLA together with the manufacturer of the drug.

Fine Mesh Baskets

We are also able to manufacture fine mesh baskets. Baskets up to 100 Mesh can be manufactured simply by replacing the mesh. For higher (finer) baskets, such as 150 mesh, the screen is too fragile to be used on its own and a supporting 20 mesh is used on the outside. This process reduces the chances of accidental damage and extends the working lifetime of the basket. Baskets are available for any bath; contact this office for a quotation.

Kioldóásvizsgálati szűrőkről, további tartozékokról részletes információ:
<http://www.labhut.com/products/tablet-dissolution.html>